



Technical Data Sheet

Streptavidin-RPE Conjugate

Package

Streptavidin-RPE is supplied in 0.1 M Sodium-phosphate buffer, pH7.2, containing 0.09% sodium azide.

Application

Streptavidin-RPE can be excited at 488 nm (absorbance max. 565 nm) with an emission maximum at approximate 575 nm.

Streptavidin-RPE conjugate could be used as an avidin/biotin labeling for Flow Cytometry, FIA and Immuno-histo staining.

Instruction

Mixed well before use. For maximal recovery of contents, please quick-spin vial before opening.

Since variety of applications, we suggest you conduct dilution test for best result determination.

Diluted sample only prepared prior use. For research or further manufacturing use only.

Storage

Store Streptavidin-RPE in dark at 2~8°C . Do not freeze.